

## Before Use

- Biosafety cabinet on?
- Drain valve closed?
- All waste containers and supplies placed inside the biosafety cabinet?
- Disinfectant prepared and in biosafety cabinet?
- Interior of biosafety cabinet disinfected?

## During Use

- Perform all work inside the biosafety cabinet
- Keep front and rear grills clear
- Discard waste inside the biosafety cabinet
- Disinfect all items prior to removal
- Remove gloves or, if wearing nitrile gloves, disinfect them before removing hands from the biosafety cabinet

## Work Completion

- Seal all waste containers before removal from the biosafety cabinet
- Disinfect all items before removal
- Disinfect all work surfaces upon completion: back, sides, inside front view screen, rear grill, work surface and front grill